




Reparaturzeichnung • Repair services drawings • Eclatés pour les services de réparation • Esplosi per servizi d'assistenza
Diseños para servicios de reparaciones • Загрузка чертежей Для ремонтных мастерских

1440-xxxx Waxprofi

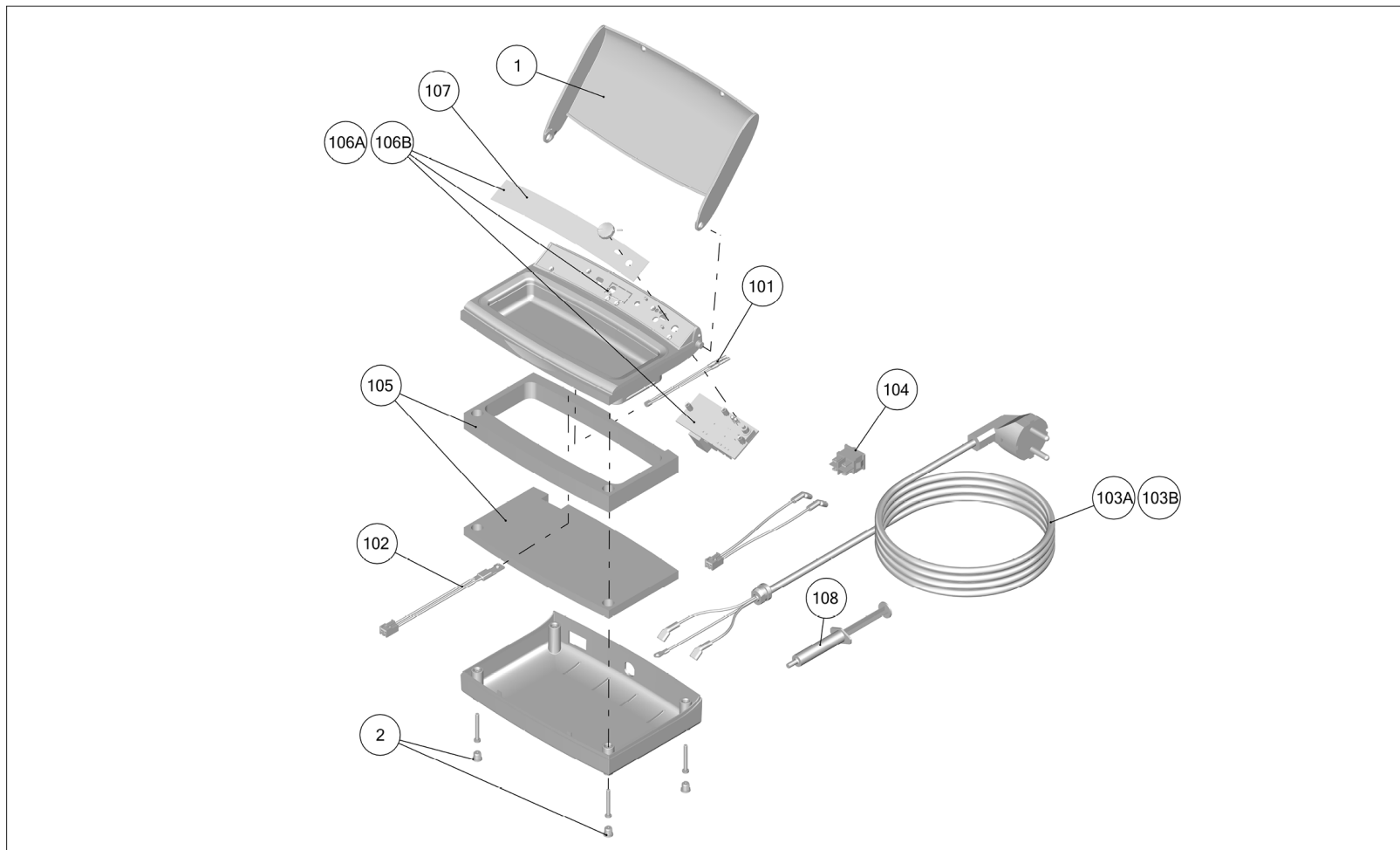

21-9903 05062018 / A



Version	Position	Name	Article number	Wishlist
A	1	Lid RAL 7035	900034543	
A	2	Rubber bases	900034564	
A	101	Thermo sensing device complete	900035783	
A	102	Thermofuse 1440	900020486	
A	103A	Power cord (230 V)	900035762	
A	103B	Power cord (120 V)	900035763	
A	104	Single-throw double-pole switch	900035747	
A	105	Isolation Set Waxprofi 1440	900034573	
A	106A	Printed boards 1440 (230V)	900034576	
A	106B	Printed boards 1440 (120V)	900034577	
A	107	Front label Waxprofi 1440	900034575	
A	108	Thermoconducting paste siliconfree	900031955	

*) Wear and tear parts, excluded from the guarantee.